



Aerodyne Mini-TILDAS CO₂ Isotope Monitor for $\Delta^{17}\text{O} - \text{CO}_2$

**Direct Spectroscopic Measurement
of $\Delta^{17}\text{O} - \text{CO}_2$ with No Chemical
Processing or Separation.**




Features:

- < 0.15 ‰ precision for $\delta^{18}\text{O}$ in 1 s.
- < 0.15 ‰ precision for $\delta^{17}\text{O}$ in 1 s.
- Fast time response (10 Hz).
- Direct measurement of CO₂ isotopes in air without sample processing.
- Repeatability exceeding 0.03 ‰ for $\Delta^{17}\text{O}$ for a 30-minute measurement including balanced working reference measurements.
- Suitable for CO₂ samples derived from carbonate via acid digestion.

TILDAS TECHNOLOGY

Aerodyne instruments use tunable infrared laser direct absorption spectroscopy (TILDAS) at mid-IR wavelengths to probe molecules at their strongest “finger-print” transition frequencies. We further enhance sensitivity by employing a patented multi-pass broad-band absorption cell that provides optical path lengths up to 76 m. Direct absorption spectroscopy allows for fast (<1 sec) absolute trace gas concentrations without need for elaborate calibration procedures. Moreover, TILDAS instruments are relatively free of measurement interference from other molecular species, enabling extremely specific detection.

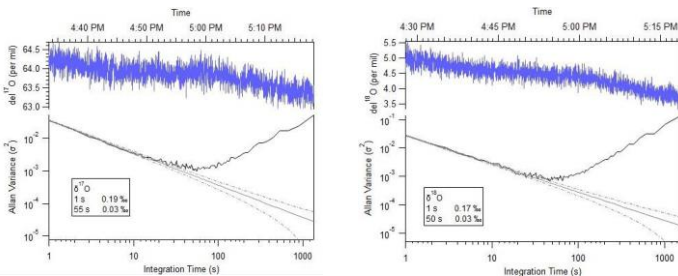

Instrument precision for a column density equivalent to 400ppm CO₂ mixing ratio and 36 meter optical path length.

Rugged, field-ready instruments

Direct absorption spectroscopy allows for highly specific and accurate gas detection

Mid-IR detection enables maximum measurement sensitivity

APPLICATIONS

- Determination of atmospheric sources, sinks, and transport through CO₂ isotopic ratios.
- Biosphere exchange.
- Analysis of CO₂ samples derived from marine carbonate samples.
- Stratosphere-troposphere exchange.

AERODYNE CO₂ ISOTOPE ADVANTAGES

- Direct measurement of ¹⁷O-CO₂ which is not possible by IRMS.
- Optional automated sample handling systems.
- Time response up to 10 Hz enables eddy covariance studies.
- Powerful TDLWintel software provides flexible instrument control, and real-time data analysis.
- Valve control capable of complex scheduling, and automatic backgrounds and calibrations.
- 19” rack mountable for easy installation.
- Turn-key design allows unattended operation in remote field sites.

CO₂ Isotope Monitor for $\Delta^{17}\text{O} - \text{CO}_2$

SPECIFICATIONS

Continuous Air Measurement Specifications – High Speed

	CO ₂	$\delta\text{CO}^{18}\text{O}$	$\delta\text{CO}^{17}\text{O}$
0.1 Second	0.15 ppm	0.6 ‰	0.6 ‰
1 Second	0.05 ppm	0.2 ‰	0.2 ‰
60 Second	0.015 ppm	0.04 ‰	0.04 ‰

Note: These measurements are not referenced to a working reference. This configuration supports 10 Hz eddy covariance measurements with a modest sized pump (120 lpm) and a flow rate of 6 slpm.

Discrete Air Sample Specifications – High Precision

	CO ₂	$\delta\text{CO}^{18}\text{O}$	$\delta\text{CO}^{17}\text{O}$
One air sample (~15 ml). 3 min measurement	0.03 ppm	0.06‰	0.06‰
10 Air samples (total of ~150ml) 30 min measurement	0.01 ppm	0.02‰	0.02‰

Dynamic Range (air)

	min	max
CO₂	0 ppm	1,000 ppm

Related Instruments

Single laser isotope monitor for $\delta^{13}\text{C}$ and $\delta^{18}\text{O}$ of CO₂

Dual laser isotope monitor for $\delta^{13}\text{C}$, $\delta^{18}\text{O}$ and $\Delta^{17}\text{O}$ of CO₂

Dual laser monitor for CO₂ ($\delta^{13}\text{C}$, $\delta^{18}\text{O}$) and water ($\delta^{18}\text{O}$, δD) isotopes

Aerodyne specializes in collaboration and custom design. Please contact us if you would like to discuss additional measurement options and applications.

REFERENCES

Wehr, R., Munger, J.W., McManus, J.B., Nelson, D.D., Zahniser, M.S., Davidson, E.A., Wofsy, S.C. and Saleska, S.R., Seasonality of temperate forest photosynthesis and daytime respiration. *Nature*, 534 (7609), pp.680-683, 2016.

Sakai, S., Matsuda, S., Hikida, T., Shimono, A., McManus, J.B., Zahniser, M., Nelson, D., Dettman, D.L., Yang, D. and Ohkouchi, N., High-Precision Simultaneous $^{18}\text{O}/^{16}\text{O}$, $^{13}\text{C}/^{12}\text{C}$, and $^{17}\text{O}/^{16}\text{O}$ Analyses for Microgram Quantities of CaCO₃ by Tunable Infrared Laser Absorption Spectroscopy. *Analytical chemistry*, 89(21), pp.11846-11852, 2017.

McManus, J. Barry, David D. Nelson, and Mark S. Zahniser. Design and performance of a dual-laser instrument for multiple isotopologues of carbon dioxide and water, *Optics express* 23.5, 6569-6586, 2015.

Wehr, R., et al. Long-term eddy covariance measurements of the isotopic composition of the ecosystem-atmosphere exchange of CO₂ in a temperate forest." *Agricultural and forest meteorology* 181, 2013, 69-84, 2013.

Installation

19" rack mountable or benchtop
Flushing the optics with CO₂-free gas is recommended

Instrument Operations

Operating temperature: 10 to 35 °C
Sample flow rate: 0 to 20 slpm

Instrument Components

Core instrument
Thermoelectric chiller
Keyboard, mouse, and monitor
Vacuum pump (customer specified)
Inlet sampling system (customizable)

Data Outputs

RS-232, USB drive, ethernet

Size, Weight, Power

Dimensions: 440 mm x 660 mm x 6U (267mm) (W x D x H)
Weight: 35 kg (core instrument) + 15 kg (chiller) + pump weight
Electrical Power: 250 W, 120/240 V, 50/60 Hz (without pump)